CL SERIES

Instruction Manual

Thank you for purchasing CL SERIES pocket scale.
Please read all operating instructions carefully before use. This electronic scale is a precision instrument. With normal care and proper treatment, it will provide years of reliable service.

Manual- English

NOTE:

- *Never load the scale with more than the maximal capacity.
- *Avoiding any exposure to extreme heat or cold, your scale works better when operated at normal room temperature. Always allow the scale to acclimate to normal room temperature for at least one hour before use.
- *Give your scale sufficient warm up time. Usually 30-60 seconds before calibrating the scale will be sufficient for most applications.
- *Keep your scale in a clean environment. Dust,dirt,misture,vibration,air currents and a close distance to other electronic equipment can all cause an adverse effect on the reliability and accuracy of your scale.
- *Handle with care. Gently approach items to be weighed onto tray top. After weighing, gently remove to be sure during try to avoid rough treatment as this may permanently damage the internal sensor.
- *Avoid shaking, dropping or otherwise shocking the scale. This is a precision instrument and must be handled with extreme care.
- *Only operate the scale on a stable, vibration free surface.

1

I. OPERATION

- Place scale on a horizontal flat surface
1. Press the [U] key to turn the scale on.
2. Wait until "0.0" or "0.00" is displayed.
3. Place the object(s) on the weighing platform.
4. By pressing the [M] key, you can switch between weighing units.

II. TARE WEIGHING

1. Turn on the scale.
2. Place a tray or container on the scale platform.
3. Press [T] key and wait until "0.0" or "0.00" is displayed
4. Add the "net- weight-item".

III. CALIBRATION:

1. Turn on the scale.
2. Press and hold the [M] key until the LCD display "CAL", release the [MODE] key.
3. Press [M] key again, the display will show "CAL" followed by the required calibration weight value.
4. Place the required calibration weight on the center of the platform, the display will show "PASS" then return to the normal weighing mode.
Calibration is complete! You can turn off the scale.

IV. ERROR CODES

1. O-LD - Maximal capacity has been exceeded. Remove your load immediately to avoid damage to the load cell.
2. L - Batteries are low. Replace batteries.
3. OUTZ - The zero range has shifted on your scale. Calibrate your scale as instructed in Section III. If the problem occurs after proper calibration, the load cell may be permanently damaged as a result of overloading.

V. SPECIFICATIONS

Tare range: Tare full capacity
Operation temperature: 10-30°C

Capacity: 300g/4629.6gn/1500oz/192.9dwt/10.583oz/9.645ozt
Readability: 0.01g/0.15gn/0.005oz/0.01dwt/0.001oz/0.001ozt
Required calibration weight: 100g/200g.
Units: g/oz/ozt/oz/ozt/gt

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